ASM-I Media Preparation Directions

The media used in this laboratory (Sasner and Haney) is based upon the ASM-I media from; Charmichael, W. and Gorham, P.; 1974, Journal Of Phycology 10: 238-240. This media is based upon a media from; Gorham, Maclachlan, Hammer and Kim,; 1964, Ver. Int. Ver. Theor. Angu. Limn. 15: 798-804 (there should be a copy of both these papers in this notebook). The chemical constituents and their concentrations for each of these media are

The chemical constituents and their concentrations for each of these media are as follows in mg/l.

Chemical	Gorham 1964	Gorham74	Sasner/Haney
NaNO ₃	170.2	170	169.8
MgCl2*6H2O	40.7	19.2	100.8
MgSO4*7H2O	49.3	24.08	86.3
CaCl ₂ *2H ₂ O	22.2	22.2	29.0
K ₂ HPO ₄	17.42	17.4	17.4
Na ₂ HPO ₄ *7H ₂ O	14.2	14.2	50.63
FeCl ₂ *6H ₂ O	1.88	.65	1.8
MnCl2*4H2O	3.89	.87	2.175
ZnCl ₂	.436	.44	1.63
CoCl2*6H2O	.0019	.01	.00348
CuCl2*2H2O	.00013	6 .0001	.000172
H ₂ BO ₃	2.47	2.47	2.47
Na ₂ EDTA*2H ₂ O		6.64	8.24

There are some variations to the media used in the Sasner/Haney cultures. the predominate ones being an increase in the amounts of magnesium, sodium/phosphate and Zinc. I could find no documentation as to the reason for this and have assumed that a less alligitrophic media was needed. In any case this is the recipe used.

The first 4 chemicals are combined into one stock media solution to from a Stock Minerals Solution. To get a 5ml stock solution for 1 liter of media you put the following amounts of chemicals into your stock solution.

Chemical	Sasner/Haney	(g/1)
NaNO ₃	33.96	
MgCl ₂ *6H ₂ O	20.16	
MgSO4*7H2O	17.36	
CaCl2*2H2O	5.8	

The next 2 chemicals are combined to form the Stock Phosphate Solution. For a 5ml stock solution into 1 liter media use the following amounts.

Chemical	Sasner/Haney	(g/1)
K ₂ HPO ₄	3.58	
Na ₂ HPO ₄ *7H ₂ O	10.12	

The next five chemicals are combined to form the Stock Metals Solution. For a 5ml stock solution into 1 liter media use the following amounts.

Chemical	Sasner/Haney	
	·	
FeCl ₂ *6H ₂ O	.36	
MnCl2*4H2O	.43	
ZnCl ₂	.33	
CoCl2*6H2O	.0069	
CuCl ₂ *2H ₂ O	.00034	
H ₂ BO ₃	.49	

The last chemical is made up separately and used as a buffer in the final media to bring it to a pH of approximately 7 to 7.5.